[탁상용 연마기 & 광택기 Grinder & Polisher BGR-10-2]

■ Characteristics

- For grinding and polishing works of rock thin plate
- Cooling water pump for internal circulation
- Possible of speed control dependent on the hardness of rock
- Aluminum plate body
- Grinder plate of stainless steel (for rust prevention)
- Magnetic plate attachment
- Grinding by diamond disc (roughly 10 times higher grinding speed than abrasive material)

■ Specification

Model		BGR - 10 - 2
Size	mm	800(W)x670(D)x230(H)
Motor		1/4HP, 220V, 2P
Plate	mm	250*2
Diamond Disc	mm	250
RPM		0-400
Cooling Pump		220V, 7W, 1.7m 800ℓ/hr

■ Accessories

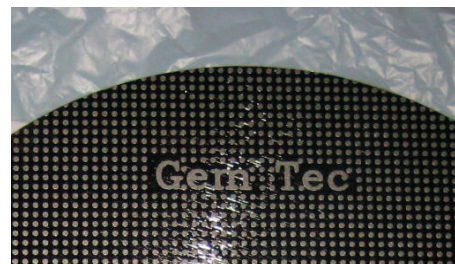
- Cooling Water Pump: 1 set
- Dot Type Diamond Disc: 200# 1 ea
- Optional disc: 2000, 400, 600, 800, 1500, 2000#
- Optional Slide Glass Holder: 28x48mm <The below photo>

<Dot Type Diamond Disc>

<Enlargement of the Topside>

<Merits of Dot Type, As Compared with the Conventional Type>

- Faster polishing speed
- Clean working environment due to the rapid releasing of the grinded ones
- Better durability
- <Patent pending>

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This equipment is a device for finely grinding the rock piece before adhering it to the slide glass.

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This equipment grinds rock piece using diamond disks or diamond cloth.

[탁상용 연마기 & 광택기 Grinder & Polisher BGR-10-2]

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Preparations before using equipment

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Check if the coolant pump contains coolant (cooling water + coolant oil) (fill 2/3 of the coolant tank).

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The coolant pump is installed at the bottom of the box.

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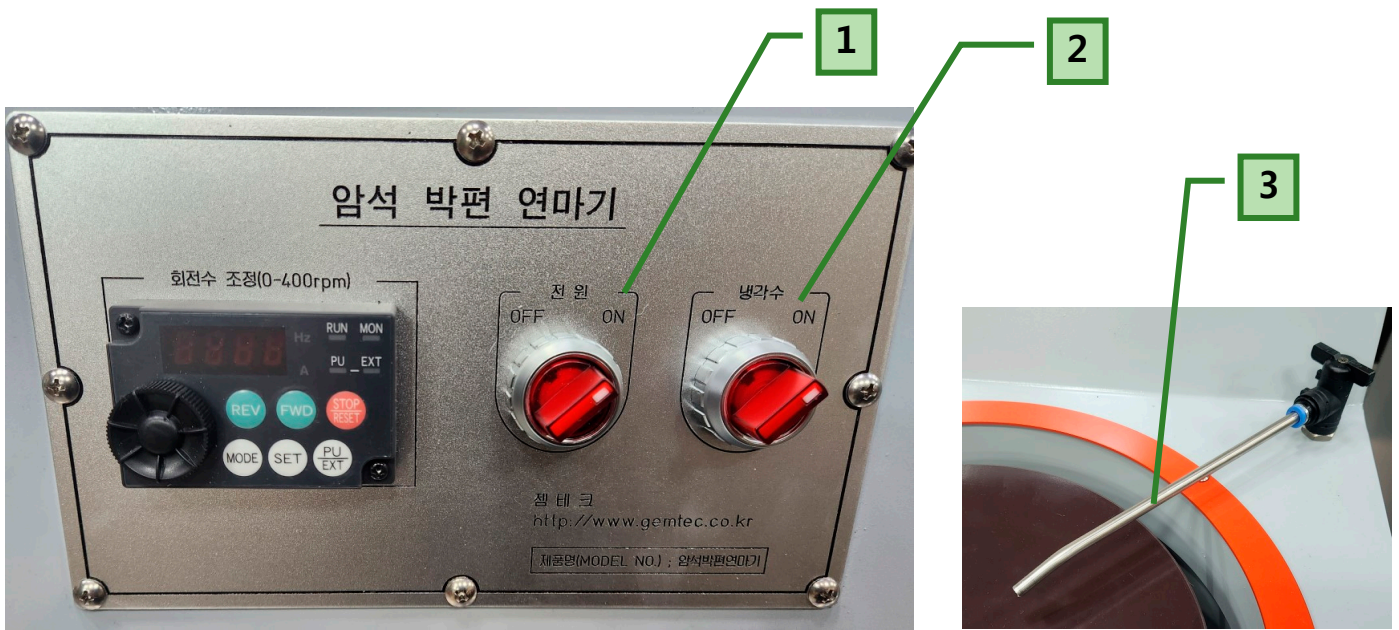
Mix the water-soluble coolant oil with water (the ratio is water 30 to oil 1, 30 vs 1). If there is no oil, you can use clean water.)

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Adjust the valve hose to supply coolant to the disc.

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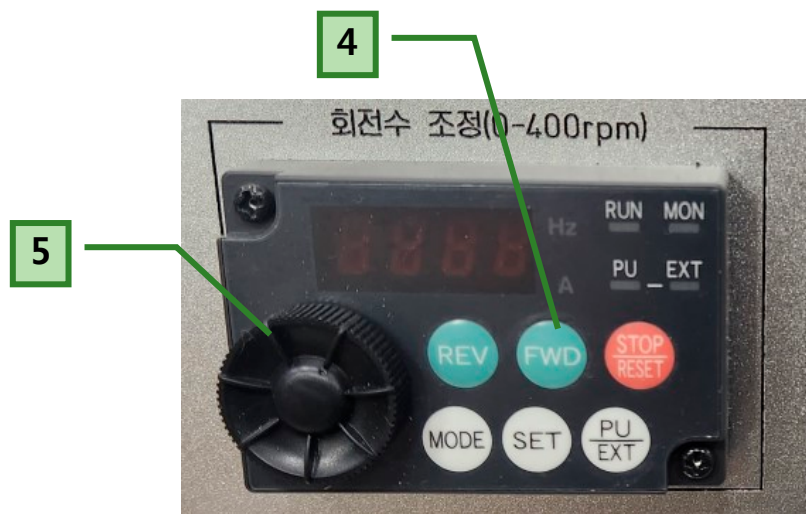
Adjust the cutting speed by adjusting the circular lever left and right. (Rotation adjustment according to specimen's strength).

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Discard the coolant after use and fill it with new coolant the next time the equipment is used.

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Use a diamond disc for the magnetic plate.

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Use an aluminum plate for silicon carbide (SiC) paper.

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When replacing the aluminum plate, turn the plate to the right to lock it and turn it to the left to unlock it. Be sure to turn off the power before replacement.

[탁상용 연마기 & 광택기 Grinder & Polisher BGR-10-2]

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Safety device for emergency stop during operation

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If you press the emergency stop (**비상정지**) button during operation, the power is cut off and all functions are turned off.

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In order for operation again, turn the emergency stop (**비상정지**) button to make the button come out.

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Switch on the switches of 1-Power, 2-Coolant.

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