[폴리셔 Polisher BGR-10]

■ Characteristics

- For grinding and polishing works of rock thin plate
- Cooling water pump for internal circulation
- Possible of speed control dependent on the hardness of rock
- Aluminum plate body
- Grinder plate of stainless steel (for rust prevention)
- Magnetic plate attachment
- Grinding by diamond disc (roughly 10 times higher grinding speed than abrasive material)

■ Specification

Model		BGR - 10
Size	mm	550(W)x550(D)x600(H)
Motor		1/4HP, 220V, 2P (380V,3P)
Plate	mm	205
Diamond Disc	mm	205
RPM		0-800
Cooling Pump		220V,7W 1.7m 800l/hr

■ Accessories

- Cooling Water Pump: 1 set
- Dot Type Diamond Disc: 200# 1 ea
- Optional disc: 2000, 400, 600, 800, 1500, 2000#
- Optional Slide Glass Holder: 28x48mm <The below photo>

<Dot Type Diamond Disc>

<Enlargement of the Topside>

<Merits of Dot Type, As Compared with the Conventional Type>

- Faster polishing speed
- Clean working environment due to the rapid releasing of the grinded ones
- Better durability
- <Patent pending>

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This equipment is a device for finely grinding the rock piece before adhering it to the slide glass.

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This equipment grinds rock piece using diamond disks or silicon paper.

[폴리셔 Polisher BGR-10]

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Preparations before using equipment

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Check if the coolant pump contains coolant (cooling water + coolant oil) (fill 2/3 of the coolant tank).

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The coolant pump is installed at the bottom of the box.

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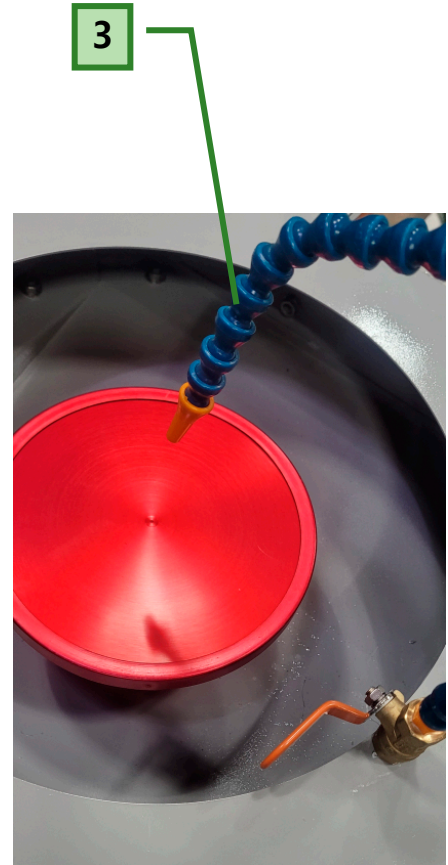
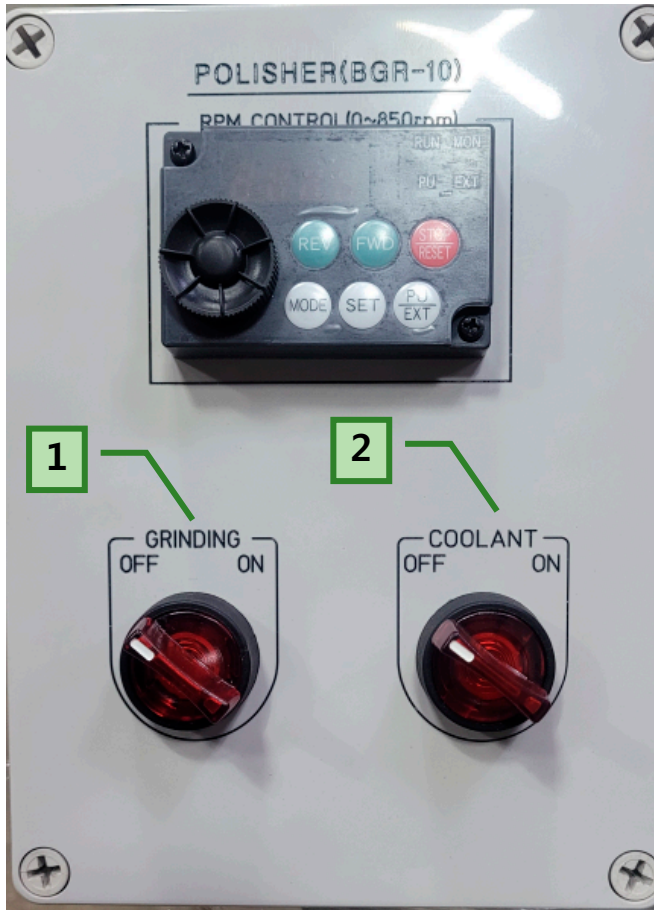
Mix the water-soluble coolant oil with water (the ratio is water 30 to oil 1, 30 vs 1). If there is no oil, you can use clean water.)

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Adjust the cutting speed by adjusting the circular lever left and right. (Rotation adjustment according to specimen's strength).

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Discard the coolant after use and fill it with new coolant the next time the equipment is used.

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Use a diamond disc for the magnetic plate.

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Use an aluminum plate for silicon carbide (SiC) paper.

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When replacing the aluminum plate, turn the plate to the right to lock it and turn it to the left to unlock it. Be sure to turn off the power before replacement.