## [암석분쇄기 Ring Mill RRM-10]

### ■ Characteristics

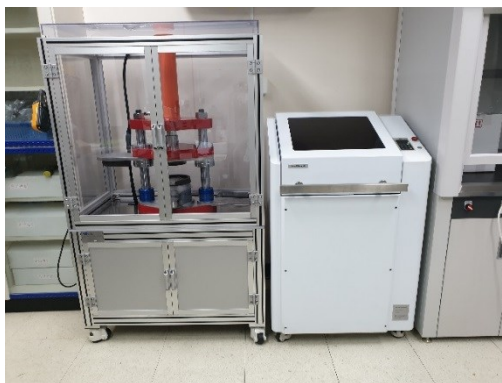
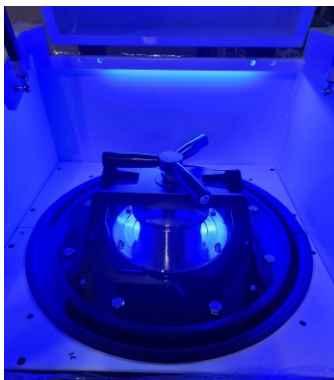
- For making rock or mineral powder with regular mesh size
- Automatic stop after powdering work
- Timer for powdering time adjustment
- Adjustment of vibration strength for powdering
- LED lighting for powdering work observation during operation
- Use of tungsten carbide to protect rock powder from any contamination
- Device for attachment and de-attachment of head air
- Case for noise-proof
- Open type air cylinder

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This equipment is a device for finely grinding the rock or mineral sample to a regular mesh size.

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This equipment uses agate, zirconia, or tungsten carbide to protect the rock or mineral sample from contamination.

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**Target group:** Operator, All persons concerned with the machine in any form.

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The RRM-10 ring mill is used for the fast, loss-free and reproducible pulverization of medium hard, hard, brittle and fibrous materials to analytical fineness. The ring mill runs smoothly and safely, also with heavy grinding sets and at maximum speed.

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The ring mill is used successfully in almost all areas of industry and research.

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This applies in particular where there are high demands in terms of hygiene, speed, fineness and reproducibility.

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The RRM-10 is particularly suitable for preparing samples for spectral analysis because it achieves high degrees of final fineness in the shortest time.

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## [암석분쇄기 Ring Mill RRM-10]

Due to its robust design, the RRM-10 has proven to be especially good in the construction materials sector (cement), in geology, mineralogy, metallurgy and in power plants.

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Soil, concrete, electronic components, ores, glass, ceramics, coal, coke, corundum, metal oxide, minerals, plant parts, slag, silicates, cement, cement clinker and many other substances can be ground easily, fast and without loss.

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This machine is a laboratory machine designed for 8-hour single-shift operation.

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This machine may not be used as a production machine nor is it intended for continuous operation.

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Increased wear or damage to the grinding set is possible if the grinding set is operated without filling or with insufficient filling quantity.

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**The grinding set must be filled to at least 1/3 of the nominal volume.**

◆ The above product may be modified for upgrade without any prior notices. ◆

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In addition to the instrument settings, the filling level of the grinding jar is also of crucial importance for a successful grinding process in the Ring Mill. The table indicates approximate values for the recommended amount of sample as a function of the grinding set used. A median filling (sample quantity) results in best grinding efficiency and increases the life time of the grinding sets.

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The mains connection must be fused to at least 16A.

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The presence of abrasive composite materials during grinding can considerably reduce tool service life.

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**When grinding electronic scrap, take the properties of the composite materials into account.**

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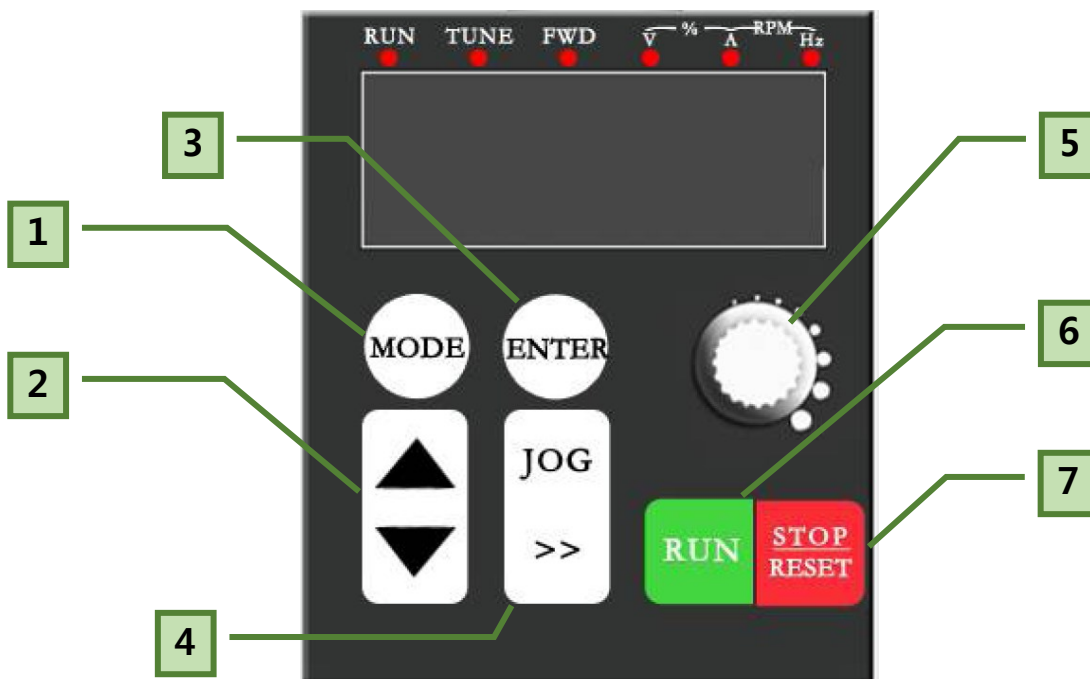
- 1500min<sup>-1</sup>

The speed can be adjusted within the following range:

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Press it to start a programs modification or cancel the modification.

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All modes can be selected and parameters adjustable with these two buttons.

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Inching run mode when the machine is in stopped state. Move the cursor to the position where needs to be modified.

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Connect the machine to the mains supply.

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Switch on the machine with the main switch at the right of the machine.

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There should be nothing in the grinding jar support.

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The VD600 can be used with a grinding set with a volume of 50 / 100 / 250ml / 500ml and with stainless steel / agate / zirconia / tungsten carbide materials.

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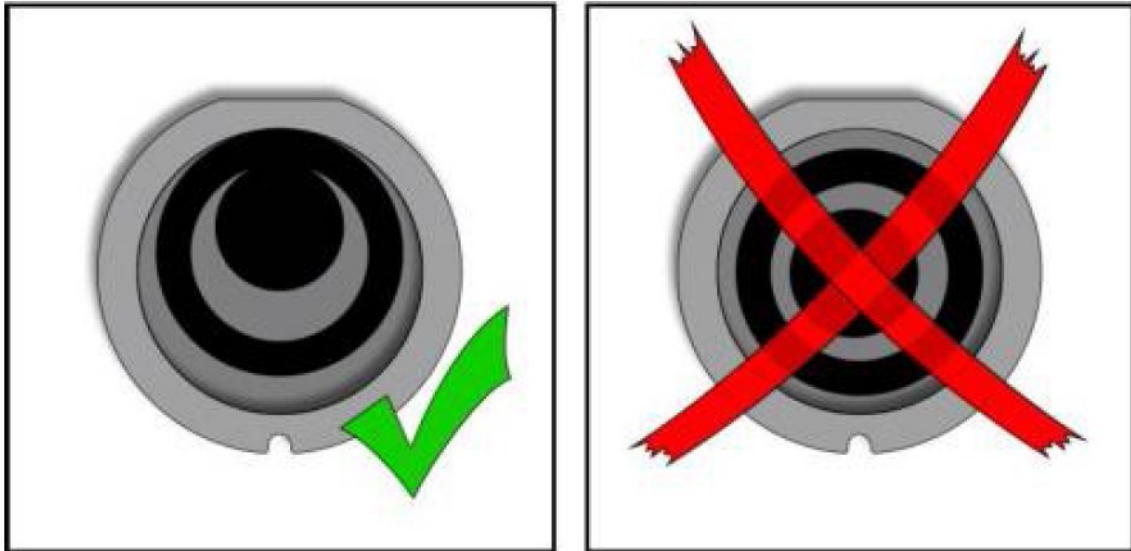
Place a disc and a ring on the grinding jar.

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Position the ring and disc as shown in the diagram.

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
Fill grinding material into the space (RZ) between the grinding jar wall (JW) and the inner ring (IR).

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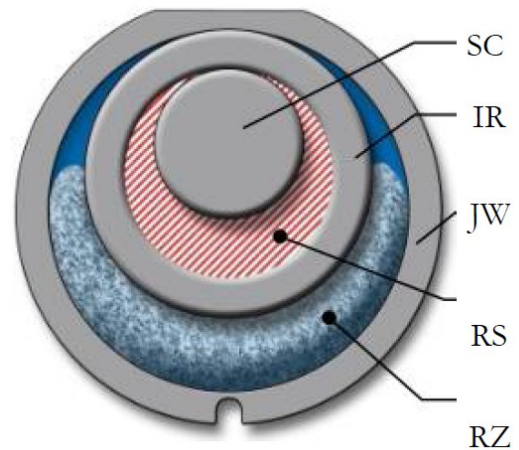For best grinding results, the area (RS) between the inner ring (IR) and the disc (SC) must remain empty.

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Remove any residual grinding material on the grinding jar wall, the inner ring and the disc, e.g. using a brush.

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Check whether the lid seal has been inserted.

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Push the grinding jar into the support.

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Clamp the grinding set using the clamping lever.

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You can then close the hood and start the device.

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Depending on the grinding process, the material being ground and accordingly the grinding jar can become very hot.

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Wear appropriate protection always when touching the grinding jar if it is hot.

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Never open hot grinding jars! Allow the grinding jars to cool down to room temperature before opening.

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The grinding chamber clamping mechanism is easy to use and reliable. The basic prerequisite both for safety of the operator and for the life of the machine components is the accurate clamping of grinding jars.

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Please remember that this device concerns grinding equipment which applies very high amounts of energy into the sample material, and that the grinding set therefore needs to be locked into position carefully.

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- The device can only be started when the clamping lever has been clamped.

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You can adjust a countdown to the starting of the device here.

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Press the STOP button to cancel the countdown.

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You can save previously set parameters such as grinding duration and speed in a memory here.

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It will show 1500, this refer to the grinding speed.

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Press [ENTER] to modify the grinding duration.

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After the modification, press [ENTER] to save the parameter.

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Finally to press [MODE] to quit program modification.

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An electric shock can cause injuries in the form of burns and cardiac arrhythmia, respiratory arrest or cardiac arrest.

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Do not clean the blender under running water. Use only a cloth dampened with water.

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Disconnect the power supply plug before cleaning the blender.

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The device must always be switched off and disconnected from the mains before any interventions for cleaning or servicing purposes

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Strong voltage through capacitor discharge

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Due to capacitor discharge on the frequency inverter, the device conducts voltage for up to 3 minutes after the plug has been pulled out.

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You may come into contact with live contacts when the device is open.

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An electric shock can lead to burns and arrhythmia or to respiratory failure and heart failure.

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