

Installation of the Operating Software for Stylus Tip Wear Observation (For the CERCHAR Abrasion Tester Microscope, Model UM22)

page: 1/2


1. Operating Software Installation

Note:

Please install the operating software on a computer or laptop equipped with USB ports and running a recent version of Windows. On older versions such as Windows 7, the software may install but may not operate properly.

Installation Steps:

- Visit the official website: <http://www.vitiny.com/en>
- On the top menu bar, open the “**Downloads**” dropdown menu.
- Locate the **UM22** model in the product list as follows:

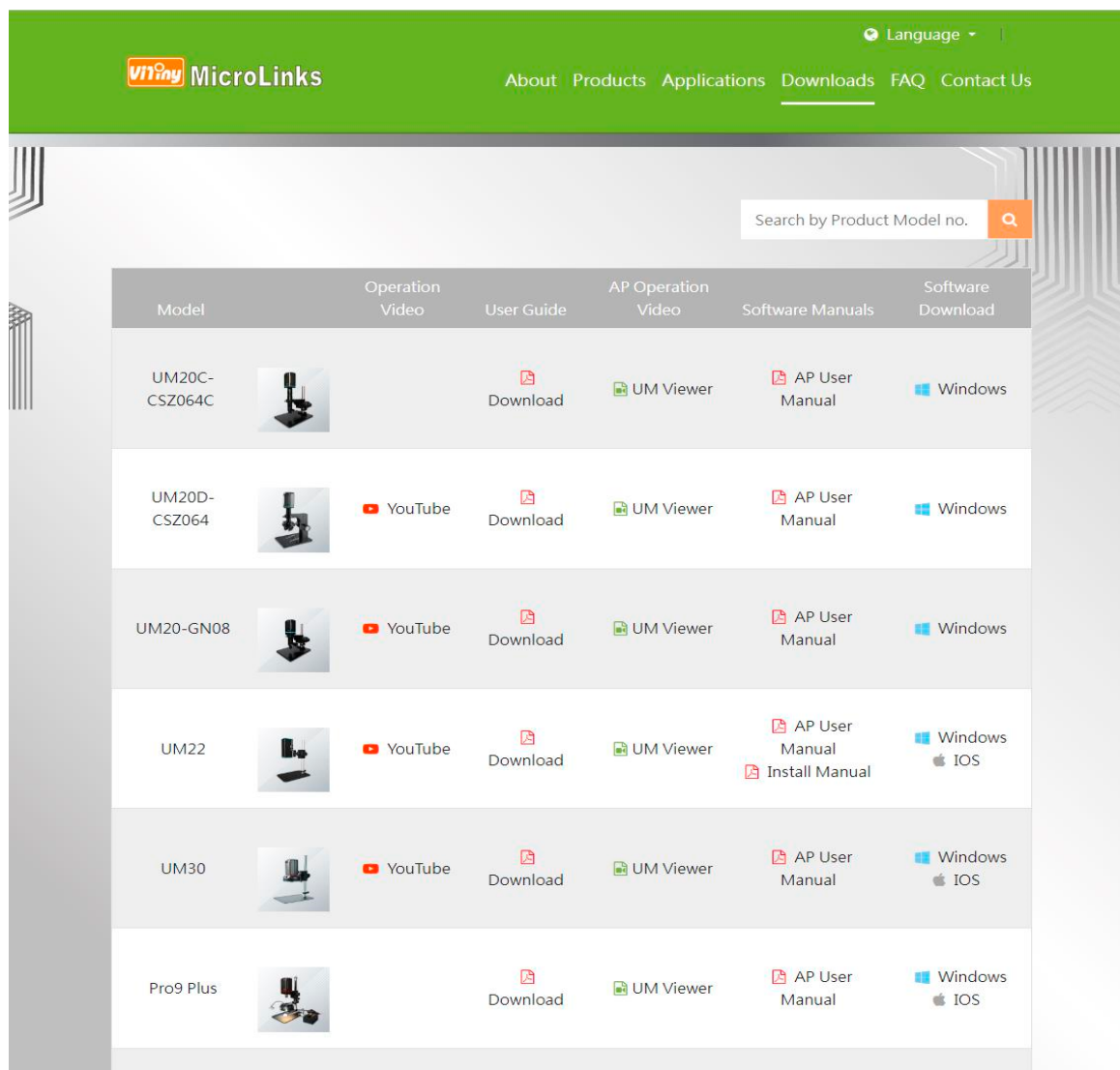

Model	Operation Video	User Guide	AP Operation Video	Software Manuals	Software Download
UM20C-CSZ064C		 Download	 UM Viewer	 AP User Manual	 Windows
UM20D-CSZ064	 YouTube	 Download	 UM Viewer	 AP User Manual	 Windows
UM20-GN08	 YouTube	 Download	 UM Viewer	 AP User Manual	 Windows
UM22	 YouTube	 Download	 UM Viewer	 AP User Manual  Install Manual	 Windows  iOS
UM30	 YouTube	 Download	 UM Viewer	 AP User Manual	 Windows  iOS
Pro9 Plus		 Download	 UM Viewer	 AP User Manual	 Windows  iOS

Installation of the Operating Software for Stylus Tip Wear Observation (For the CERCHAR Abrasion Tester Microscope, Model UM22)

page: 2/2

- d. In the “**Application Program**” column on the far right, click “**Software Download**” to download and install the operating software.

2. Microscope-Related Documents

(Refer to the illustration above.)

- a. **User Guide:** Operating manual for the microscope
b. **Software Manuals:** Operating manual for the software

- End -